

Delta Tips

NMDT_0079


Delta V6.0 - New Save Functions

NMR data processing software





In Delta V6.0, it is possible to save multiple data displayed in the **Data Slate** window to one file in the JDA format. It is also possible to save the processed 2D data altogether with high resolution 1D data to one file in the **2D Viewer** window.


In the **1D Processor** and **Data Slate** windows :

Click the **Save All Data**  button to save all data in the **1D processor** or **Data Slate** window to a single JDA file.

★ The FID data is kept under version number 1 and the FT-processed data is stored under version numbers 2, 3, and so on.

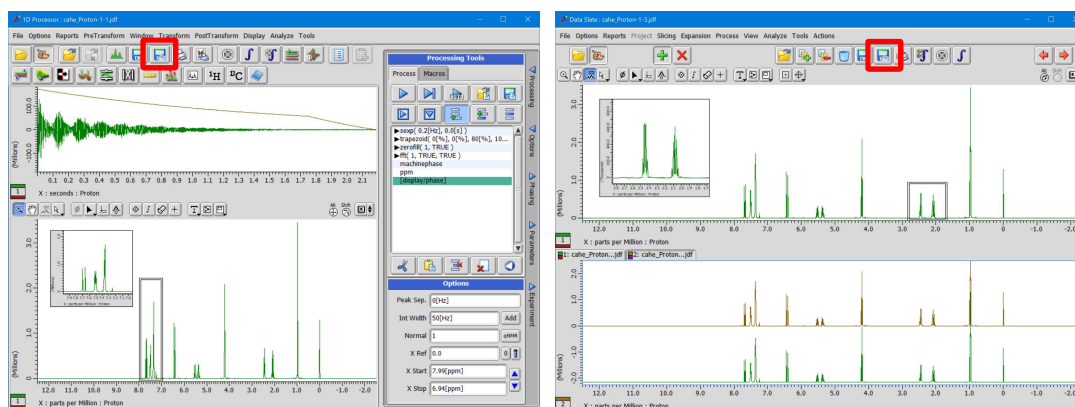


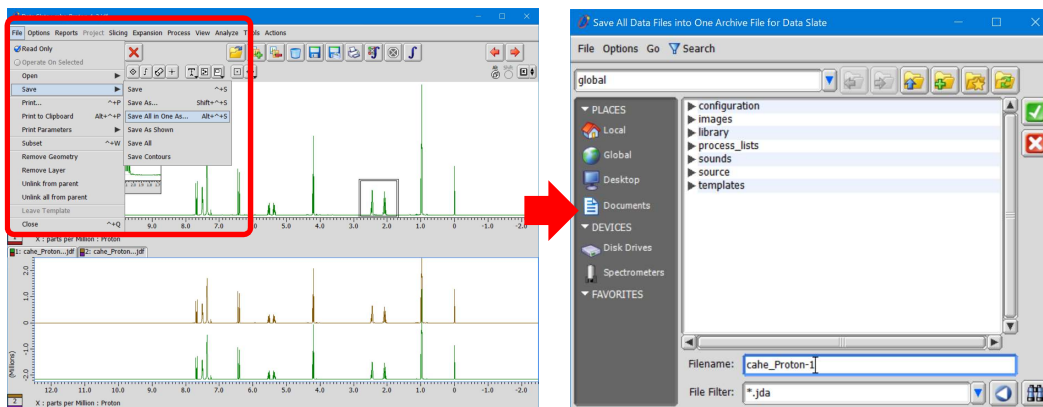
★ Click the **Save All Data**  button while holding the **Alt** key to select the directory and filename.

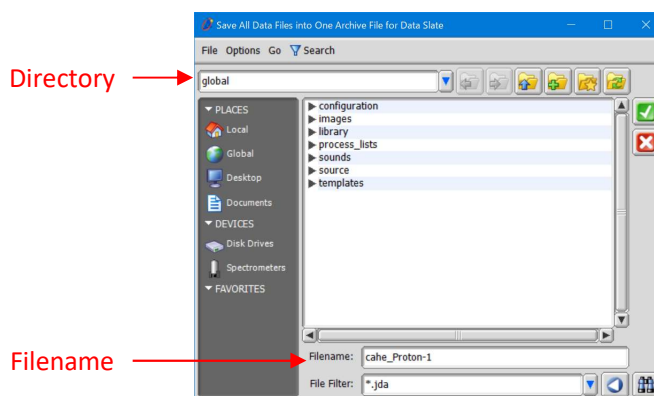

Delta Tips


- Select **File—Save—Save All in One As** in the menu bar.



This function is the same as holding the **Alt** key and clicking the **Save All Data**  button.



In the **2D Viewer** window :

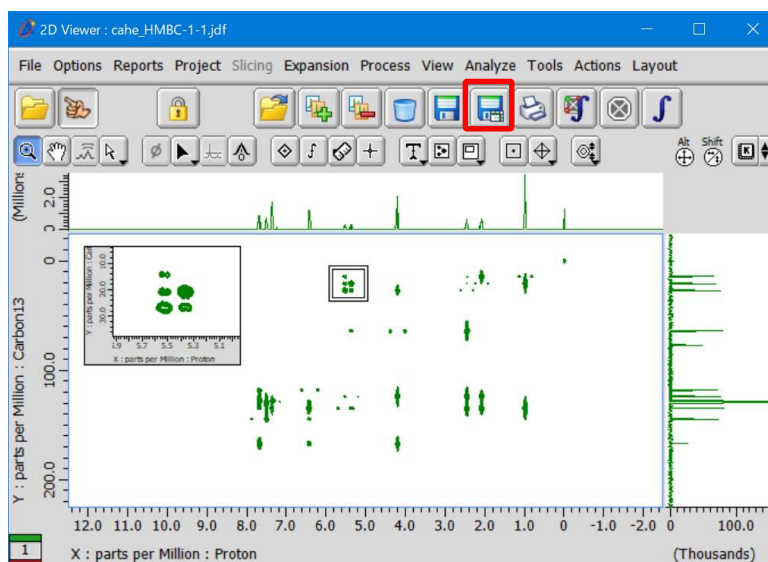Click the **Save All Data**  button to save all data altogether to a single JDA file.



- ★ In the **2D Viewer** window, the **Save All Data**  button has the same function as the **Save All Data**  button in the **Data Slate** window. It is also possible to use the combination of **Alt** key and the **Save All Data** button or select **File—Save—Save All in One as** in the menu bar.